

## Zombie NIR™ Fixable Viability Kit

<b>Catalog# / Size</b>	423105 / 100 tests 423106 / 500 tests
<b>Regulatory Status</b>	RUO
<b>Other Names</b>	Fixable Dye, Fixable Viability Dye, Dead, Apoptosis
<b>Description</b>	Zombie NIR™ is an amine reactive fluorescent dye that is non-permeant to live cells, but permeant to the cells with compromised membranes. Thus, it can be used to assess live vs. dead status of mammalian cells. Zombie NIR™ is a polar water soluble dye, providing red fluorescence, making it suitable for multi-color detection.

### Product Details

---

<b>Preparation</b>	Zombie NIR™ Fixable Viability kit is composed of lyophilized Zombie NIR™ dye and anhydrous DMSO. For reconstitution, bring the kit to room temperature; add 100 µl of DMSO to one vial of Zombie NIR™ dye until fully dissolved. 100 tests = 1 vial of Zombie NIR™ + DMSO, 500 tests = 5 vials of Zombie NIR™ + DMSO.
<b>Storage &amp; Handling</b>	Store kit at -20°C upon receipt. Do not open vials until needed. Once the DMSO is added to the Zombie NIR™ dye, use immediately, or store at -20°C in a dry place and protected from light, preferably in a desiccator or in a container with desiccant for no more than one month.
<b>Application</b>	<a href="#">FC - Quality tested</a>
<b>Recommended Usage</b>	Each lot of this product is quality control tested by immunofluorescent staining with flow cytometric analysis.  For flow cytometry, the suggested dilution is 1:100-1:1000 for 1-10 million cells. It is recommended that the reagent be titrated for optimal performance for each application, as optimal dosage varies with cell type.
<b>Excitation Laser</b>	Red Laser (633 nm)
<b>Application Notes</b>	Zombie NIR™ dye is excited by the red laser and has fluorescence emission maximum at 746 nm. If using in a multi-color panel design, filter optimization may be required depending on other fluorophores used. Zombie NIR™ dye has similar emission to APC/Cyanine7.

#### Standard Cell Staining Protocol:

1. Prior to reconstitution, spin down the vial of lyophilized reagent in a microcentrifuge to ensure the reagent is at the bottom of the vial.
2. For reconstitution, pre-warm the kit to room temperature; add 100 µL of DMSO to one vial of Zombie NIR™ dye and mix until fully dissolved
3. Wash cells with PBS buffer (no Tris buffer and protein free).
4. Dilute Zombie NIR™ dye at 1:100-1000 in PBS. Resuspend 1-10 x 10<sup>6</sup> cells in diluted 100 µL Zombie NIR™ solution. To minimize background staining of live cells, titrate the amount of dye and/or number of cells per 100 µL for optimal performance. Different cell types can have a wide degree of variability in staining based on cell size and degree of cell death.
  1. **Note:** Don't use Tris buffer as a diluent and be sure that the PBS does not contain any other protein like BSA or FBS.
  2. **Note:** The amount of dye used can also influence the ability to detect apoptotic as well as live and dead cells.
5. Incubate the cells at room temperature (or 4°C), in the dark, for 15-30 minutes.
6. Wash one time with 2 mL BioLegend's Cell Staining Buffer (Cat. No. 420201) or equivalent buffer containing serum or BSA.
7. Continue performing antibody staining procedure as desired.

8. Cells can be fixed with paraformaldehyde or methanol prior to permeabilization or can be analyzed without fixation.

#### No-wash Sequential Staining Protocol:

1. Wash cells with PBS buffer (no Tris buffer and protein free).
2. For reconstitution, pre-warm the kit to room temperature; add 100  $\mu$ L of DMSO to one vial of Zombie NIR™ dye and mix until fully dissolved
3. Determine the total  $\mu$ L volume of antibody cocktail previously titrated and optimized for the assay that will be added to each vial/well of cells based on a final volume of 100  $\mu$ L. Subtract that antibody volume from the 100  $\mu$ L total staining volume intended for the assay. In the remaining volume, dilute Zombie NIR™ dye at 1:100-1000 in PBS as determined by prior optimization at that volume. For example, if you are adding 20  $\mu$ L of antibody cocktail for a 100  $\mu$ L total staining volume, use 80  $\mu$ L of Zombie NIR™ solution. Resuspend  $1-10 \times 10^6$  cells in the appropriate volume of Zombie NIR™ solution. Different cell types can have a wide degree of variability in staining based on cell size and degree of cell death.

1. **Note:** Don't use Tris buffer as a diluent and be sure that the PBS does not contain any other protein like BSA or FBS.
2. **Note:** The amount of dye used can also influence the ability to detect apoptotic as well as live and dead cells.

4. Incubate for 10-15 minutes at RT (or 4°C), protected from light. Without washing the cells, add the cell surface antibody cocktail and incubate for another 15-20 minutes.
5. Add 1-2 mL Cell Staining Buffer (Cat. No. 420201) or equivalent buffer containing BSA or serum. Centrifuge to pellet.
6. Continue with normal fixation and permeabilization procedure. If planning to skip fixation and analyze cells live, complete an additional wash step to minimize any unnecessary background of the live cells.

1. **Notes:** If the cell type in use cannot tolerate a protein-free environment, then titrate the Zombie NIR™ dye in the presence of the same amount of BSA/serum as will be present in the antibody staining procedure. A higher amount of Zombie NIR™ may be required since the BSA/serum will react with and bind up some proportion of the Zombie NIR™.

**Additional Product Notes** View more applications data for this product in our [Scientific Poster Library](#).

#### Application References

(PubMed link indicates BioLegend citation)

1. McMaster SR, *et al.* 2015. *PLoS One*. 10:115725. [PubMed](#)
2. Prado-Garcia H, *et al.* 2015. *Anticancer Res*.35:1529. [PubMed](#)
3. Rodriguez-Rodriguez N, *et al.* 2015. *J Immunol*. 194:4207. [PubMed](#)
4. Flies DB, *et al.* 2015. *J Immunol*. 194:5294. [PubMed](#)

#### Product Citations

1. Komuczki J, *et al.* 2019. *Immunity*. 50:1289. [PubMed](#)
2. Rad S MAH, *et al.* 2020. *PLoS One*. 15:e0232915. [PubMed](#)
3. Headley M, *et al.* 2016. *Nature*. 531:513-517. [PubMed](#)
4. Wang L, *et al.* 2020. *Front Cell Dev Biol*. 0.745138889. [PubMed](#)
5. Iyer S, *et al.* 2021. *Cancer Discov*. 0.725. [PubMed](#)
6. Kim MY, *et al.* 2021. *JCI Insight*. 6:. [PubMed](#)
7. Martínez-Vélez N, *et al.* 2022. *Mol Cancer Ther*. 21:471. [PubMed](#)
8. Baratchart E, *et al.* 2022. *PLoS Comput Biol*. 18:e1009839. [PubMed](#)
9. Luo J, *et al.* 2022. *J Nanobiotechnology*. 20:228. [PubMed](#)
10. Rosina M, *et al.* 2022. *Cell Metab*. 34:533. [PubMed](#)
11. Hsu MA, *et al.* 2022. *Cancer Immunol Immunother*. :. [PubMed](#)
12. Zheng J, *et al.* 2022. *Front Pharmacol*. 13:866228. [PubMed](#)

## Antigen Details


---

**Biology Area** Apoptosis/Tumor Suppressors/Cell Death, Cell Biology, Neuroscience

**Gene ID** NA

## Product Data

---

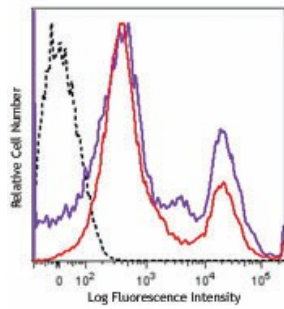

One day old C57BL/6 mouse splenocytes were stained with Zombie NIR™ and analyzed before fixation (purple) or after fixation and permeabilization (red). Cells alone, without Zombie NIR™ staining, are indicated in black.

For Research Use Only. Not for diagnostic or therapeutic use.

This product is supplied subject to the terms and conditions, including the limited license, located at [www.biolegend.com/terms](http://www.biolegend.com/terms) ("Terms") and may be used only as provided in the Terms. Without limiting the foregoing, BioLegend products may not be used for any Commercial Purpose as defined in the Terms, resold in any form, used in manufacturing, or reverse engineered, sequenced, or otherwise studied or used to learn its design or composition without express written approval of BioLegend. Regardless of the information given in this document, user is solely responsible for determining any license requirements necessary for user's intended use and assumes all risk and liability arising from use of the product. BioLegend is not responsible for patent infringement or any other risks or liabilities whatsoever resulting from the use of its products.

BioLegend, the BioLegend logo, and all other trademarks are property of BioLegend, Inc. or their respective owners, and all rights are reserved.

8999 BioLegend Way, San Diego, CA 92121 [www.biolegend.com](http://www.biolegend.com)  
Toll-Free Phone: 1-877-Bio-Legend (246-5343) Phone: (858) 768-5800 Fax: (877) 455-9587